



Sterile, alcohol-based ready to use solution for disinfecting surfaces in cleanrooms.

## perform<sup>®</sup> sterile alcohol EP

### Our Plus

- fast and effective
- sterile (γ-irradiated) and microbially filtered (0.2 μm)
- suitable for cleanrooms grade A/B, ISO class 5/6
- ready to use in a handy, sterile 500 ml spray system
- colour and perfume free
- efficacy tested according to European norms (EN)

### Application areas

Disinfection of surfaces and equipment/devices in cleanrooms. Especially for Grade A and B clean areas, according to:

- EU Guide to Good Manufacturing Practice for medicinal products Annex 1: Manufacture of sterile medicinal products, March 2009.
- FDA Draft Guidance for Industry: Sterile Drug Products produced by aseptic processing - current Good Manufacturing practice, U.S. Department of Health and Human Services; Food and Drug Administration, September 2004.

### Instructions for use

Apply perform<sup>®</sup> sterile alcohol EP undiluted to objects and surfaces and allow to take effect. Ensure complete wetting. The product dries off quickly. perform<sup>®</sup> sterile alcohol EP can be used for glove disinfection. Maximum application rate 50 ml/sq.m. Do not use to treat surfaces that are particularly sensitive to alcohol, such as acrylic glass.

### Microbiological efficacy

Efficacy	Concentration	Contact time
bactericidal EN1276, EN13697	ready-to-use	1 min.

Efficacy	Concentration	Contact time
fungicidal EN1650, EN13697	ready-to-use	2 min.
virucidal against enveloped viruses in accordance with DVV (German Registered Association for Combating Viral Diseases)//RKI Guideline	ready-to-use	30 sec.



# perform<sup>®</sup> sterile alcohol EP

## Product data

Composition:

100 g solution contains the following active substances: 25 g Ethanol (94 %), 35 g Propan-1-ol.

## Special advice

**Use biocides safely. Always read the label and product information before use.**

For all of the related safety information please refer to the Safety Data Sheet. To request the Safety Data Sheet please E-mail: [tech.uk@schuelke.com](mailto:tech.uk@schuelke.com)

The sterile 500 ml spray system offers the following advantages:

- convenient size for storage
- ergonomically shaped spray pump for ease of use
- remains sterile after opening

IE: PCS Number 95504

## Information for order

Item	Delivery form	Item no.
perform sterile alcohol EP 500 ml FL	10/Carton	121802

## Environmental information

schülke manufactures products economically and with advanced, safe and environmentally friendly production processes while at the same time maintaining out high quality standards.

## Expert opinion and information

Please visit our website for an overview of all available literature/reports on all schulke products online at <http://www.schuelke.com/>

For any queries you might have about the product or using it please call 0114 2543500 or E-mail [mail.uk@schuelke.com](mailto:mail.uk@schuelke.com)



Schülke & Mayr GmbH holds a Manufacturer's Authorisation according to sect 13 para 1 German Drug Law and Certificates of GMP Compliance for medicinal products.



**schülke UK**  
Schülke & Mayr UK Ltd  
Cygnet House, Meadowhall  
Sheffield, S9 1AT  
Phone +44-114-254 35 00  
Fax +44-114-254 35 01  
[mail.uk@schuelke.com](mailto:mail.uk@schuelke.com)

**schülke Headquarters**  
Schülke & Mayr GmbH  
Robert-Koch-Str. 2  
22851 Norderstedt, Germany  
Phone +49 40 - 52100 - 0  
Fax +49 40 - 52100 - 318  
[www.schuelke.com](http://www.schuelke.com)  
[info@schuelke.com](mailto:info@schuelke.com)