



Microbially filtered, alcohol-based ready to use solution for disinfecting surfaces in cleanrooms.

perform® advanced alcohol EP

Our Plus

- fast, effective and residue free
- microbially filtered (0.2 µm)
- ready to use
- suitable for cleanrooms grade C/D
- colour and perfume free
- efficacy tested according to European norms (EN)

Application areas

Disinfection of surfaces and equipment/devices in clean room areas and other areas with strict hygienic requirements.

Instructions for use

Apply perform® advanced alcohol EP undiluted to objects and surfaces and allow to take effect. Ensure complete wetting. The product dries off quickly. perform® advanced alcohol EP is also suitable for disinfecting gloves. Maximum application rate 50 ml/sq.m. Do not use to treat surfaces that are particularly sensitive to alcohol, such as acrylic glass.

Microbiological efficacy

Efficacy	Concentration	Contact time
bactericidal <i>EN1276, EN13697</i>	ready-to-use	1 min.
fungicidal <i>EN1650, EN13697</i>	ready-to-use	2 min.
virucidal against enveloped viruses <i>in accordance with DVV (German Registered Association for Combating Viral Diseases)//RKI Guideline</i>	ready-to-use	30 sec.

Efficacy	Concentration	Contact time
Vaccinia virus <i>in accordance with DVV (German Registered Association for Combating Viral Diseases)//RKI Guideline</i>	ready-to-use	30 sec.



perform[®] advanced alcohol EP

Product data

Composition:

100 g solution contains the following active substances: 25 g Ethanol (94 %), 35 g Propan-1-ol.

Special advice

Use biocides safely. Always read the label and product information before use.

For all of the related safety information please refer to the Safety Data Sheet. To request the Safety Data Sheet please

E-mail: tech.uk@schuelke.com

IE: PCS Number 95497

Information for order

Item	Delivery form	Item no.
perform advanced alcohol EP 1 l FL	10/Carton	121804
perform advanced alcohol EP 10 l KA	1/Canister	121805

Environmental information

schulke manufactures products economically and with advanced, safe and environmentally friendly production processes while at the same time maintaining out high quality standards.

Expert opinion and information

Please visit our website for an overview of all available literature/reports on all schulke products online at <http://www.schuelke.com/>

For any queries you might have about the product or using it please call 0114 2543500 or E-mail mail.uk@schuelke.com



Schülke & Mayr GmbH holds a Manufacturer's Authorisation according to sect 13 para 1 German Drug Law and Certificates of GMP Compliance for medicinal products.

schulke UK
Schülke & Mayr UK Ltd
Cygnet House, Meadowhall
Sheffield, S9 1AT
Phone +44-114-254 35 00
Fax +44-114-254 35 01
mail.uk@schuelke.com

schulke Headquarters
Schülke & Mayr GmbH
Robert-Koch-Str. 2
22851 Norderstedt, Germany
Phone +49 40 - 52100 - 0
Fax +49 40 - 52100 - 318
www.schuelke.com
info@schuelke.com