



Sterile, ready to use 70 % (v/v) IPA impregnated wipes for surface disinfection in cleanrooms.

perform® sterile wipes IPA

Our Plus

- fast, effective and residue free
- pouch of 20 ready to use alcohol impregnated wipes
- sterile (X-Ray irradiated) and double wrapped
- suitable for cleanrooms grade A/B, ISO class 5/6
- low particle release
- colour and perfume free
- efficacy tested according to European norms (EN)

Application areas

Disinfection of surfaces and equipment / devices in cleanrooms, in particular for class A and B cleanroom areas, according to:

- Annex 1 of EU Guide to Good Manufacturing Practice Manufacture of sterile medicinal products, March 2009.
- FDA Guidance for Industry: Sterile drug products produced by aseptic processing - current Good Manufacturing Practice; U.S. Department of Health and Human Services, Food and Drug Administration, September 2004.

Instructions for use

For professional use only. Clean the surface carefully before use. Remove excess water from the surface before disinfection, if necessary. Ensure complete wetting. Used wipes must be disposed of in a sealed container. Disinfect after each production and cleaning process or if required according to SOP. Wipe the surface thoroughly with a cloth and leave at room temperature (20± 2°C) for at least 5 minutes. Application for PT02: Disinfection of clean non-porous surfaces such as small work surfaces in medical and non medical areas as well as surfaces in clean rooms (class A/B). Effective against bacteria and yeasts.

Microbiological efficacy

Efficacy	Concentration	Contact time
bactericidal EN1276 - clean conditions	ready-to-use	5 min.
bactericidal EN16615 - clean conditions	ready-to-use	5 min.
yeastocidal EN1650 - clean conditions	ready-to-use	5 min.
yeastocidal EN16615 - clean conditions	ready-to-use	5 min.



perform[®] sterile wipes IPA

Product data

Composition:

100 g solution contains the following active substances: 63.1 g Propan-2-ol (equates 70% (v/v)).

Chemical-physical data

Color	colourless
Density	ca. 0,88 g/cm ³ / 20 °C
Flash point	22 °C / Method : DIN 51755 Part 1
Form	liquid
pH	Not applicable
Viscosity, dynamic	> 2,5 mPa*s / 20 °C / Raw material

Special advice

Use disinfectants safely. Always read the label and product information before use.

The alcohol used within the perform[®] sterile wipes IPA comprises alcohol of pharmaceutical quality in accordance with the Ph. Eur.

EU-0023656-0002 1-2

Information for order

Item	Delivery form	Item no.
perform [®] sterile wipes IPA 20 pc bag	20/Carton	on request

These products are not available in every country. For more information please contact our local subsidiary or distributor.

Related Products

- perform[®] sterile alcohol EP
- perform[®] sterile alcohol IPA
- perform[®] sterile alcohol IPA (WFI)
- perform[®] sterile dry wipes
- perform[®] sterile mix dry wipes
- perform[®] sterile premium dry wipes

Environmental information

schülke manufactures products economically and with advanced, safe and environmentally friendly production processes while at the same time maintaining our high quality standards.

Expert opinion and information

Please visit our website for an overview of all available literature/reports on the product: www.schuelke.com. For individual questions:

Customer Sales Service

Phone: +49 40 52100-666

E-Mail: info@schuelke.com



Schülke & Mayr GmbH holds a Manufacturer's Authorisation according to sect 13 para 1 German Drug Law and Certificates of GMP Compliance for medicinal products.

schülke Headquarters
Schülke & Mayr GmbH
Robert-Koch-Str. 2
22851 Norderstedt, Germany
Phone +49 40 - 52100 - 0
Fax +49 40 - 52100 - 318
www.schuelke.com
info@schuelke.com