



non-sterile cellulose-polyester dry wipes suitable for use in cleanrooms

perform[®] classic mix dry wipes

Our Plus

- Low particle release
- Suitable for cleanrooms grade C/D/E
- Tested compatibility with products of the perform[®] range
- 150 wipes are single bagged

Application areas

For disinfection and cleaning activities that require the use of wipes with very low particle release. Especially suitable for use in cleanrooms grade C/D, ISO class 7/8.

Instructions for use

Apply disinfectant/cleaner on the wipe, then wipe objects and surfaces and allow to take effect. Ensure complete wetting. The compatibility of the perform[®] classic mix dry wipes was tested with all disinfectants of the perform[®] range. Special attention was given to the chemical stability of the active ingredients of the disinfectants in the presence of the wipe material.

The following products of the perform[®] range are compatible:

- perform[®] advanced alcohol EP
- perform[®] classic alcohol EP
- perform[®] classic alcohol IPA
- perform[®] classic concentrate PAA
- perform[®] classic concentrate OXY

For detailed information on test methods and parameters (e.g. contact times) please contact us.



perform[®] classic mix dry wipes

Product data

Material:

- Nonwoven hydroentangled 55% cellulose and 45% polyester wipe

Size:

- 30 x 30 cm

Thickness:

- 0,27 - 0,34 mm

Packaging:

- 150 wipes are single bagged. One carton contains 10 of such bags with 150 wipes.

Surface mass:

- 68 g/m²

Particle release:

- LPC ($\geq 0.5 \mu\text{m}$): $< 12 \times 10^7$ particle/m² (IEST-RP-CC 004.3 - orbital shake method)
- APC (Helmke Drum Test @10rpm); ($\geq 0.5 \mu\text{m}$): < 340 particle per wipe per m³ per hour (IES-RP-CC003.2)

Absorption (IEST-RP-CC 004.3):

- capacity: 285 ml/m²
- speed (method: drop disappearance): 0.8 seconds

Non-volatile residues (IEST-RP-CC 004.3, short extraction)*:

- Extraction with DIW (100 ml DI water, room temperature for 24 hours) < 0.2015 mg/g

Ion content (IEST-RP-CC 004.3, short extraction)*:

- Sodium (Na) < 0.04 mg/g
- Potassium (K) < 0.008 mg/g
- Chloride (Cl) < 0.05 mg/g

**(typical test results, these are not part of the product specification)*

Information for order

Item	Delivery form	Item no.
perform [®] classic mix dry wipes 150 pieces / bag	10/Carton	on request

These products are not available in every country. For more information please contact our local subsidiary or distributor.

Environmental information

schülke manufactures products economically and with advanced, safe and environmentally friendly production processes while at the same time maintaining out high quality standards.

Expert opinion and information

Please visit our website for an overview of all available literature/reports on the product: www.schuelke.com.

For individual questions:

Application Department

Phone: +49 40 52100-666

E-Mail: info@schuelke.com



Schülke & Mayr GmbH holds a Manufacturer's Authorisation according to sect 13 para 1 German Drug Law and Certificates of GMP Compliance for medicinal products.

schülke Headquarters
Schülke & Mayr GmbH
Robert-Koch-Str. 2
22851 Norderstedt, Germany
Phone +49 40 - 52100 - 0
Fax +49 40 - 52100 - 318
www.schuelke.com
info@schuelke.com